testoval[®]

Mangan - Test 0 - 20 mg/l

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Manganese Test 0 - 20 mg/l

Directions for Use

Remove the test cell from the comparator and fill it up to the mark with the water sample. Add 6 drops of reagent A and a dosing spoon of reagent B and dissolve by stirring with the red spoon.

After 1 minute, insert the test cell into the comparator. Hold the comparator against a light source or a bright background and match the colour of the central field (test solution) with one of the colours of the peripheral colour fields. Read the value shown under the matching colour field.

The colour comparison should be carried out immediately after a 1 minute waiting time. A further delay can result in the test solution turning brown, which will invalidate the test.

Should the colour in the central field be more intensive than the highest colour value in the peripheral fields, dilute the water sample 1 : 10. For this purpose, fill the graduated plastic tube up to the 1 ml mark with the water sample and dilute with manganese-free water (e.g. distilled water) to the 10 ml mark. Transfer the diluted sample into the test cell which has been rinsed previously with manganese-free water and proceed as described above. The value obtained must be multiplied by 10.

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